LINHA LS · SÉRIE X.000 · LS
LINE · SERIES X.000

Liofilizador de Bancada LS X.000

Benchtop
Freeze Dryer
LS X.000

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The LS Line consists of two main assemblies: the freeze-dryer cabinet, which houses the drying chamber, refrigeration system and control panel; and the vacuum pump, supplied separately and connected through a flexible hose.

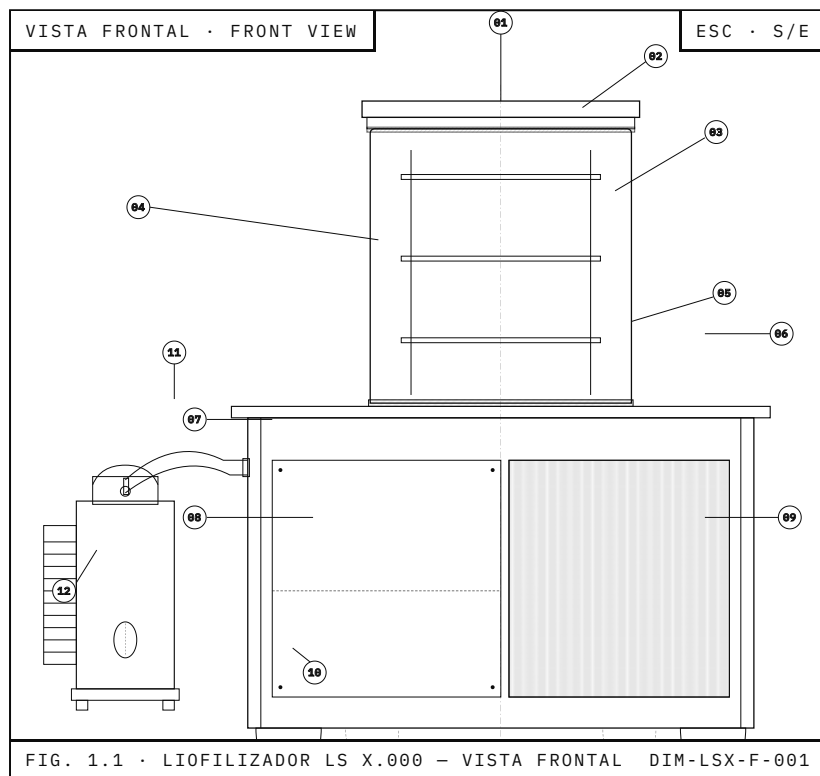
FIG. 1.1 · LIOFILIZADOR LS X.000 – VISTA FRONTAL DIM-LSX-F-001

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The LS Line is supplied in 2 capacity models — LS 3000 and LS 6000 — and 4 equipment configurations (A, B, C, D), combining tray heating and automatic vial stoppering. Check the identification label on the rear of the unit.

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D · PERFORMANCE DE PROCESSO
D · PROCESS PERFORMANCE

TEMP. CONDENSADOR CONDENSER TEMP.	-50 °C (setpoint -30 °C de fábrica) <i>(factory setpoint -30 °C)</i>
VÁCUO FINAL TÍPICO TYPICAL FINAL VACUUM	80 — 200 µHg
SETPOINT BANDEJA TRAY SETPOINT	-20 °C — +40 °C · somente A & B <i>A & B only</i>
ALARME VÁCUO VACUUM ALARM	> 300 µHg
ALARME TEMP. TEMP. ALARM	> 70 °C
TAXA DE SECAGEM DRYING RATE	± 0,8 mm / h · pH 7 (referência) <i>pH 7 (reference)</i>

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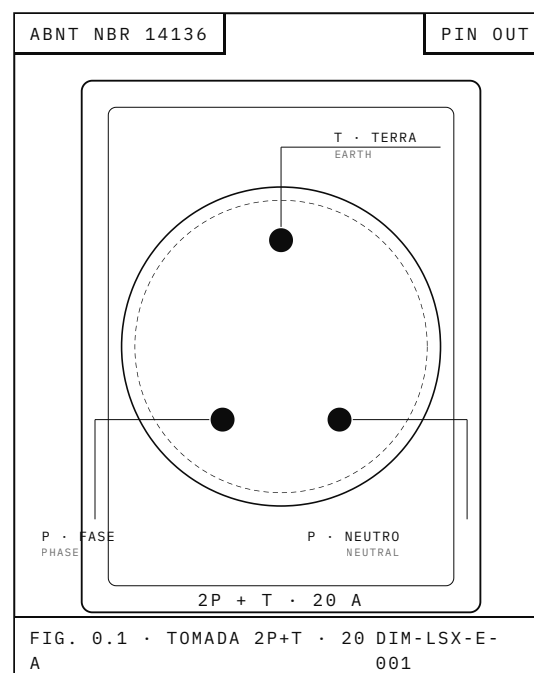
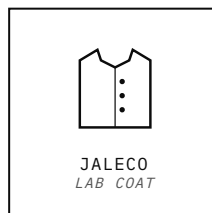
The warranty is rendered void in cases of: incorrect installation or absence of proper earthing; improper use or exposure to acid / alkaline vapours; transport damage (carrier's responsibility); connection to inadequate voltage; operation or service by personnel not authorised by the manufacturer; and natural wear of items such as seals, manifolds, acrylic parts, casters and paintwork.

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The rear panel houses every connection and identification point on the equipment. Keep it clear at least 200 mm from any wall to allow airflow through the radiator.

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Lyophilisation (freeze-drying) is the most refined process of product preservation, combining the two most reliable methods available: freezing and dehydration.

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To be lyophilised, the product must be fully crystallised (frozen through the core out to the periphery) below -20 °C . It is then exposed to negative pressure (vacuum), which removes water by sublimation — it transitions directly from solid (ice) to gas (vapour), bypassing the liquid phase.

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The final product has a porous, moisture-free structure that can be reconstituted by simply adding water. Its original properties — size, texture, flavour, aroma, vitamins, mineral salts, proteins and bacteria — are preserved and, when properly stored even at room temperature, remain intact for many years.

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If the product is not fully crystallised and is still placed in a freeze dryer, so-called vacuum dehydration occurs. Final quality is poorer, as much of the original properties are lost (they "run off" with the water).

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Total pre-freezing of the product, $\leq -20\text{ °C}$, for $\geq 24\text{ h}$ in a conventional freezer.

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Ice sublimation under vacuum. Typical rate of $\pm 0.8\text{ mm/h}$ at neutral pH.

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Removal of adsorbed moisture via controlled platform heating.

	LIOFILIZADOR LS X.000	REV 3.3	ESC S/E	LANG PT · EN	FOLIO 06 / 30
	MANUAL DE INSTRUÇÕES	DOC DIM-LSX-3.3	SEÇ. 2.0 · Intro..	APR. —	A4 297×210
DATA 05-2026	UNIT MM				

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Pre-freezing is the first and most important stage of good lyophilisation, since it directly affects the quality of the final product.

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*There is no universal rule for the freezing (crystallisation) stage, given the wide variety of products that can be lyophilised. When the product is crystallised in a conventional market freezer ($\approx -25\text{ °C}$), however, it must remain there for a **minimum of 24 hours**, provided the matrix is aqueous and the pH neutral; otherwise the time must be determined experimentally.*

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The size of crystals formed during solidification defines the appearance and many properties of the final product:

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Rapid crystallisation ($\leq -80\text{ °C}$): small ice crystals form. Dried products are lighter in colour, but small crystals tend to aggregate, hindering water-molecule migration during drying.

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Slow crystallisation: produces large ice crystals, leaving relatively open structures after sublimation — favourable for vapour migration, but large sharp crystals can rupture cellular membranes.

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Avoid freezing inside the freeze-dryer drying chamber. For most products, we recommend conventional freezers (around -20 °C) for this stage.

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The main methods to perform pre-freezing are direct-contact cooling (static) and rotational cooling in a refrigerant bath (dynamic).

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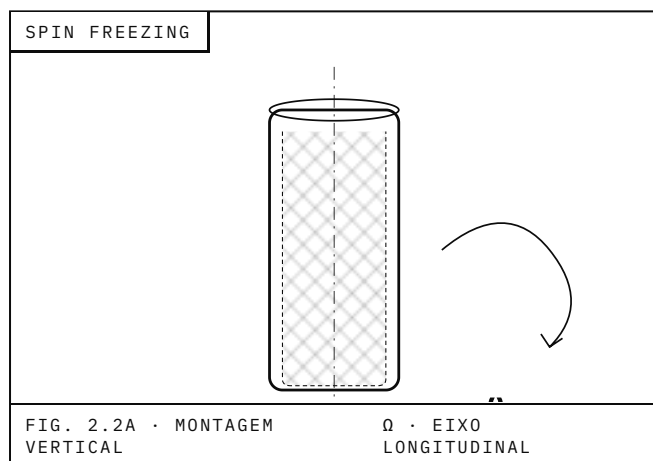
A static technique in which the cooling device drives a decrease of 1 to 4 °C/min down to the safe temperature. For most products, a conventional freezer (around -20 °C) is sufficient and recommended.

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
A dynamic method, mostly used for larger liquid volumes. It has two variants — spin freezing and shell freezing — both increase the product surface exposed to vacuum and are well-suited for vials.

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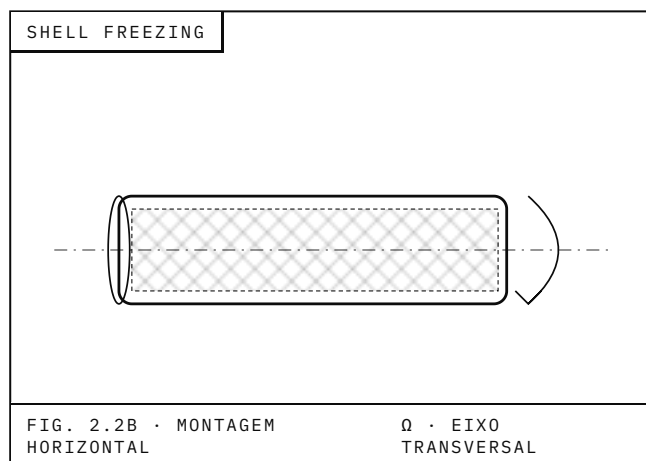The vial is mounted on a device that rotates about its longitudinal axis. Centrifugal force distributes and solidifies the product on the vial walls.

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Similar technique to spin freezing, but with horizontal mounting — the product forms a frozen “shell” on the inner walls of the container.

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Sometimes the cooled product appears fully crystallised, but it "foams" when exposed to vacuum. This happens when the core is not fully crystallised and water is pulled out by the vacuum.

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A small bubble forming on the surface of the frozen pellet indicates that the product was not fully crystallised. Despite the frozen appearance, liquid remained inside the pellet (or slab).

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The fix is to increase freezer time or reduce pellet thickness.

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Bubbles (foaming) or displacement of the pellet inside vials or trays may indicate slow transfer from freezer to drying chamber. Surfaces liquefy by air exposure during transport or by delay in starting the vacuum.

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To avoid thermal imbalance when processing small fractions, add mass by placing a water mirror in the tray or pre-cool the platforms with artificial ice packs (low-eutectic-point solutions). Remove the packs and dry the platform with paper towel before loading the product tray.

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Frozen products must look opaque and remain so throughout the drying process. After placing the product to freeze, never open the freezer to check it — if needed, wait approximately 3 hours.

	LIOFILIZADOR LS X.000	REV 3.3	ESC S/E	LANG PT · EN	FOLIO 09 / 30
	MANUAL DE INSTRUÇÕES	DOC DIM-LSX-3.3	SEÇ. 2.3 · Obser...	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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Freezing with dry ice can substitute an ultra-freezer (-40 °C to -80 °C range) for initial investment and running cost. Liquid nitrogen is recommended for products that precipitate rapidly.

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Freezing with liquid nitrogen requires special trays and extreme caution to avoid vial cracking or accidents, given its violent evaporation. Only experienced or trained personnel should perform this procedure.

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Since handling frozen vials one by one is impractical, group them on trays and move them as such from the freezer to the drying chamber. When unloading from the freezer to the chamber, work as fast as possible to prevent temperature gain and thawing during transfer.

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If only a few vials with product need to be processed, fill out the tray with empty vials, forming a thermal barrier.

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Sublimation starts at the surface of the frozen product. As the process progresses, the sublimated surface deepens into the pellet and vapours migrate through the already-dried layer. For pH 7 products the typical rate is ± 0.8 mm/h.

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Sublimation of free ice under vacuum. Most water is removed during this stage.

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Adsorbed bound water is removed. Platforms can be heated (models A and B) to speed up the process.

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For products with pH below 6 (acid) or above 8 (alkaline), the equipment must have a special vacuum pump (dry pump) or undergo frequent pump-oil decontamination/replacement.

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV 3.3	ESC S/E	LANG PT · EN	FOLIO 11 / 30
		DOC DIM-LSX-3.3	SEÇ. 2.5 · Secag..	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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The unit uses a 3-pin mains outlet (2P + earth) per ABNT NBR 14136. The centre pin must be properly earthed per the regulation below.

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Oil particle dispersion may occur through the pump exhaust. If noticeable, route the gases outdoors or to a muffler. The conduit must descend at ≥ 15° — with no siphon — and terminate outdoors (fixed PVC tube + Ø¾" translucent hose) or in a wide-mouth plastic drum containing stainless steel or copper wool.

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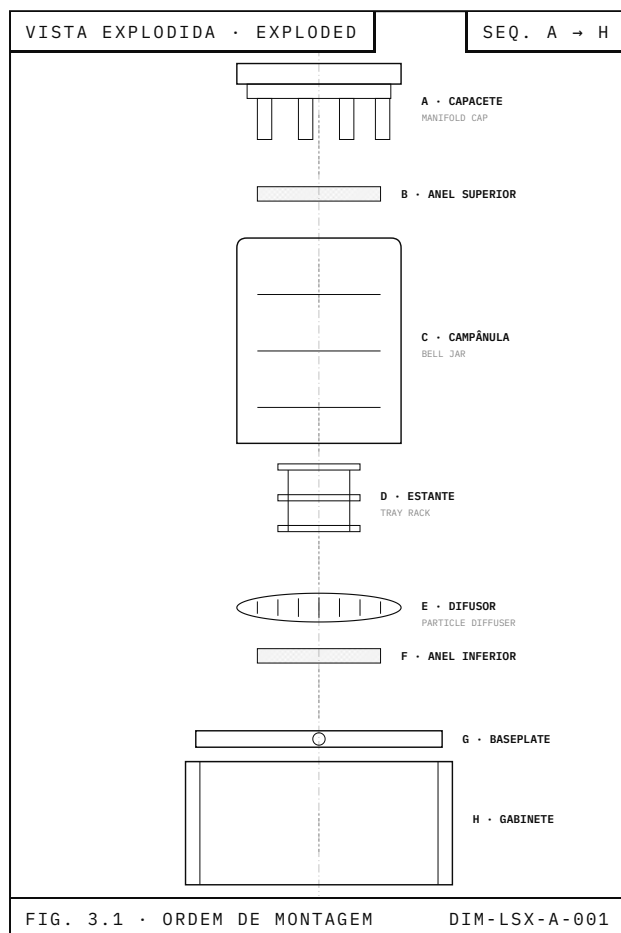
The vacuum pump is shipped separately for transport convenience. Place it beside the freeze-dryer on a sturdy bench and follow the sequence below to assemble the drying chamber.

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Always handle the bell jar with utmost care — avoid drops and damage to the acrylic structure, especially the centering on the particle diffuser.

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV 3.3	ESC S/E	LANG PT · EN	FOLIO 13 / 30
		DOC DIM-LSX-3.3	SEQ. 3.1 · Monta...	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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The unit has three menus: two external operator menus (no password) and one internal menu protected by password. On power-up the LCD shows the model and board revision; then the main status screen.

Tela principal de status *Main status screen*

VÁCUO · VACUUM	0000 µHg
CONDENSADOR · CONDENSER	00 °C
BANDEJA · TRAY	00 °C
SET POINT	00 °C
DATA · DATE	DD.MM.AA
HORA · TIME	HH:MM:SS
DECORRIDO · ELAPSED	HH:MM
REDE AC · AC LINE	000 V

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To enter: power the unit on with [01] Power and, while the version screen is visible, hold [03] ▲ until the Language parameter appears.

IDIOMA · LANGUAGE	PT · ES · EN — use [03] ▲ para alterar, [08] OK para confirmar. <i>use [03] ▲ to change, [08] OK to confirm.</i>
CONTRASTE LCD	Use [03] ▲ para aumentar, [04] ► para diminuir. <i>Use [03] ▲ to increase, [04] ► to decrease.</i>
DATA & HORA	[03] ▲ altera o valor · [08] OK confirma e avança ao próximo dígito. <i>[03] ▲ changes the value · [08] OK confirms and advances to the next digit.</i>

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV 3.3	ESC S/E	LANG PT · EN	FOLIO 15 / 30
		DOC DIM-LSX-3.3	SEÇ. 4.1 · Menus...	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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*The internal menu sets the unit's individual parameters. After powering up with [01] Power and while the version is on screen, press [02] Function — a password screen appears; enter **123 456** to access.*

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*The unit has an **RS-485** serial output and dedicated software, supplied by the manufacturer, that allows data handling, monitoring, configuration, storage, visualisation and printing through a PC.*

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Before starting, ensure the equipment is clean and dry, with no residues on the stainless steel baseplate. The same applies to the acrylic bell jar and the sealing ring.

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On models A and B with heating, the temperature sensor is on the 1st platform (bottom-up). This probe commands the heating cycle.

	LIOFILIZADOR LS X.000	REV 3.3	ESC S/E	LANG PT · EN	FOLIO 17 / 30
	MANUAL DE INSTRUÇÕES	DOC DIM-LSX-3.3	SEÇ. 5.0 · Op. C..	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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On models A and B, the system determines end-of-cycle from vacuum and platform-heating curves. Heating only starts when the platform 1 sensor reaches the start setpoint (e.g. +5 °C) and the minimum vacuum is reached (set at 200 µHg).

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*When that occurs, the system releases the heating ramp and modulates temperature per the configured power (e.g. 10%). A **lightning-bolt** icon appears on the display, indicating the platform temperature is under control.*

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At any time the user can end the cycle: press [05] Start/Stop. The screen asks "End process?" — press [08] OK to confirm or [05] to resume.

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When Tray SetP (user-defined) and Maximum Vacuum (factory: 100 µHg) are met, "Process Finished" appears.

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Without heating, consider the product ready when vacuum reaches its lowest possible level (200 — 80 µHg) and the product appears opaque, with the outside of the bell jar dry.

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Manifolds are nitrile-rubber valves designed for vial lyophilisation. They accept standard penicillin vials (Ø 20 mm) directly, or wide-mouth vials (Ø 70–80 mm) via supplied silicone adaptors.

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Before starting, check that all manifolds are in VENT position (closed to the freeze dryer) and that the condenser drain valve is closed.

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Press [05] Start/Stop to begin. The condenser cooling system engages. Upon reaching the safe temperature (-30 °C), the display reads "Place product."

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With the condenser at safe temperature, remove each vial individually from the freezer — do not connect them all at once.

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For wide-mouth vials, freeze them with the adaptor in place — handling via the silicone, never direct contact with the vial.

P05 Conecte o **bico de vidro do adaptador** à boca do manifold/válvula.

Connect the adaptor glass spout to the manifold's nozzle.

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Open the valve by turning the plunger from VENT to VAC — opens the manifold to the chamber, applying vacuum to the vial. The product must stay inert and opaque.

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Wait for vacuum recovery (back to the value before opening), then repeat for each vial until all desired manifolds are filled.

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The larger the vials used, the longer the interval between connections — and the longer to complete drying.

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV	ESC	LANG	FOLIO
		3.3	S/E	PT · EN	20 / 30
DATA	UNIT	DOC	SEÇ.	APR.	A4
05 · 2026	MM	DIM-LSX-3.3	6.0 · Manif...	—	297x210

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PROCESSAMENTO CONTÍNUO / CONTINUOUS OPERATION

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Vial processing supports continuous operation. The equipment may run all week without defrosting or draining meltwater — as long as the volume lyophilised does not exceed condenser capacity.

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*The ampoule tree connects via hose to a manifold. The remaining manifolds can run simultaneously in the same cycle. Ampoule freezing is done **externally** in a nearby freezer, and the entire tree (with ampoules attached) is then transferred.*

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV 3.3	ESC S/E	LANG PT · EN	FOLIO 21 / 30
		DOC DIM-LSX-3.3	SEÇ. 6.1 · Retir...	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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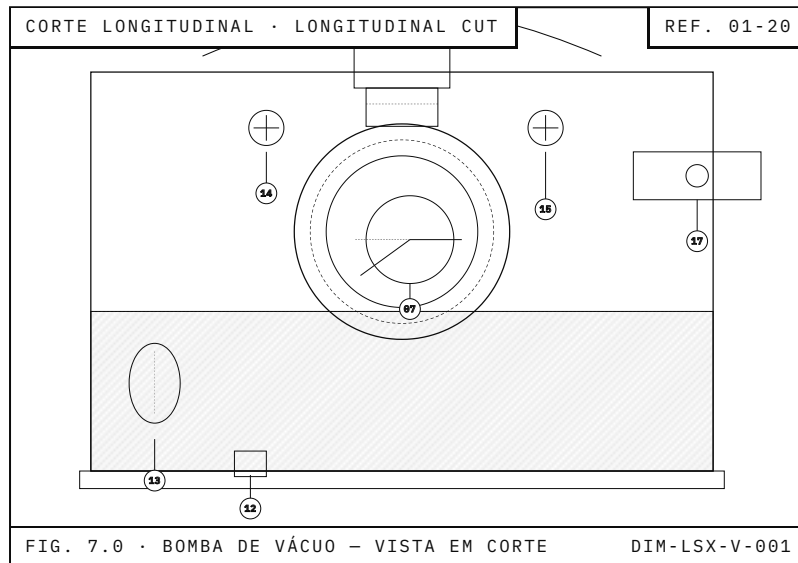
The vacuum pump is the unit's heart — oil maintenance is high priority. The better the process vacuum, the better the lyophilised sample quality. The pumping unit operates submerged in high-quality, low-vapour-pressure oil that serves as seal, coolant and lubricant.

- | | |
|----|---|
| 01 | Entrada de ar
<i>Air inlet</i> |
| 02 | Filtro
<i>Filter</i> |
| 03 | Refletor de óleo
<i>Oil deflector</i> |
| 04 | O-ring
<i>O-ring</i> |
| 05 | Mola
<i>Spring</i> |
| 06 | Palheta · vane
<i>Vane</i> |
| 07 | Rotor
<i>Rotor</i> |
| 08 | Estator
<i>Stator</i> |
| 09 | Caixa de óleo
<i>Oil box</i> |
| 10 | Óleo de vácuo
<i>Vacuum oil</i> |
| 11 | Válvula de saída
<i>Outlet valve</i> |
| 12 | Parafuso do dreno de óleo
<i>Oil drain plug</i> |
| 13 | Visor de óleo
<i>Oil sight glass</i> |
| 14 | Parafuso de entrada de óleo
<i>Oil fill plug</i> |
| 15 | Válvula de gás ballast
<i>Gas-ballast valve</i> |
| 16 | Filtro de poeira
<i>Dust filter</i> |
| 17 | Saída de gás · escapamento
<i>Gas exhaust</i> |
| 18 | Alça de transporte
<i>Carry handle</i> |
| 19 | Acoplamento flexível
<i>Flexible coupling</i> |
| 20 | Tampas de proteção
<i>Protective covers</i> |

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV 3.3	ESC S/E	LANG PT · EN	FOLIO 22 / 30
		DOC DIM-LSX-3.3	SEÇ. 7.0 · Bomba	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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Indicates sediments or acids. Immediate replacement.

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Indicates emulsified water. Immediate change.

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Indicates light contamination. Perform gas-ballast decontamination.

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With the pump running, unscrew the CAPS to open the gas-ballast valve. Air drawn through the orifice carries dissolved substances out via the exhaust, returning the oil to its original colour. Duration: 20 to 40 minutes. If the oil remains milky or dark after that period, replace it immediately. When done, close the CAPS.

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Level must always be checked with the pump running, never stopped. When stopped, oil migrates into the inner chambers and the level drops — if topped up in that state, the pump will expel the excess through the exhaust.

	LIOFILIZADOR LS X.000	REV 3.3	ESC S/E	LANG PT · EN	FOLIO 23 / 30
	MANUAL DE INSTRUÇÕES	DOC DIM-LSX-3.3	SEÇ. 7.1 · Óleo	APR. —	A4 297×210
DATA 05·2026	UNIT MM				

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The faults listed below can be addressed by the operator. Take time to understand each step and what is happening. If guidance is insufficient, contact the manufacturer's technical assistance.

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*If you encounter problems or doubts about operation, contact the manufacturer: [+XX (XX) 0000-0000] · [support@manufacturer.com]. **Do not entrust your equipment to unauthorised technicians**, even those claiming to be "specialists".*

	LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES	REV 3.3	ESC S/E	LANG PT · EN	FOLIO 24 / 30
DATA	UNIT	DOC DIM-LSX-3.3	SEÇ. 8.0 · Panes	APR. —	A4 297×210
05·2026	MM				

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The only maintenance the equipment requires is listed below. Care given to the vacuum pump is of real importance for proper operation and durability.

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Despite the externally identical look to commercial compressors, the refrigeration unit uses a special refrigerant mix — it requires critical adjustments. Do not entrust it to unauthorised technicians.

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Always keep the radiator air-intake unobstructed on the rear cabinet — clear path for circulation.

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
Care for manifold plungers: clean corrosive substances immediately — hydrochloric, sulphuric or formic acids can corrode metal parts of the chamber.

LIOFILIZADOR LS X.000 MANUAL DE INSTRUÇÕES		REV 3.3	ESC S/E	LANG PT · EN	FOLIO 25 / 30
		DOC DIM-LSX-3.3	SEQ. 9.0 · Manut...	APR. —	A4 297×210
DATA 05-2026	UNIT MM				

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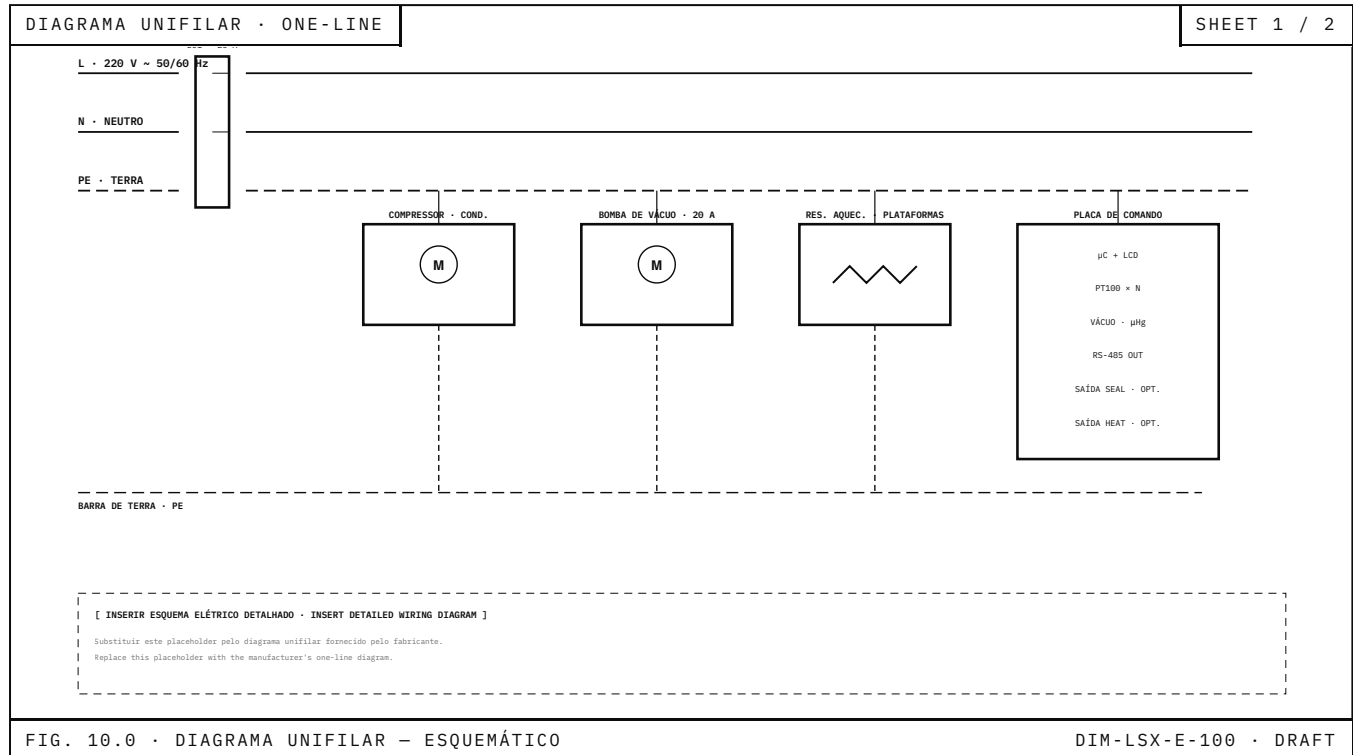Dear customer, it is with great satisfaction that we have delivered this equipment. To improve our service, please answer the questionnaire below. Your opinion is extremely important to us.

DATA ENTREGA DELIVERY DATE	
Nº PROPOSTA QUOTE NO.	
Nº NOTA FISCAL INVOICE NO.	
CLIENTE CUSTOMER	
EQUIPAMENTO · MODELO EQUIPMENT · MODEL	
RESPONSÁVEL RESPONSIBLE	

Avalie o atendimento de suas expectativas *Rate the service against your expectations*

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Avalie o atendimento de suas expectativas *Rate the service against your expectations*

	ÓTIMO GOOD	BOM GOOD	RUIM POOR
1. O técnico realizou um bom atendimento? <i>1. Did the technician deliver good service?</i>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
2. As informações foram passadas de forma clara? <i>2. Was information clearly conveyed?</i>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>
3. As dúvidas foram sanadas? <i>3. Were your questions answered?</i>	<input type="checkbox"/>	<input type="checkbox"/>	<input type="checkbox"/>

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*The information in this manual is basic guidance for understanding the process and using the equipment properly. It **may not be reproduced**, even in part, without prior written authorisation from the manufacturer. For questions, contact our Customer Service.*

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© [ANO] · [FABRICANTE] · TODOS OS DIREITOS RESERVADOS · ALL RIGHTS RESERVED